

Please refer to <http://www.medicinalgenomics.com/product-literature/> for updated protocols and Material Safety Data Sheets (MSDS). Consult MSDS before using any new product.

*PathoSEEK™ is a registered trademark of Medicinal Genomics Corporation and is for laboratory use only.*

## **Table of Contents**

PathoSEEK™ Analysis Quick Reference Tables: .....	2
FLOWER .....	2
NON – FLOWER Matrices.....	2
GUMMY .....	3
Detailed Assay Data Analysis .....	4
Table 4, Cq to CFU Conversion Equation Table.....	14
Troubleshooting Guide.....	16
Glossary and Definitions .....	17
DISCLAIMER .....	17
LIMITED USE LABEL LICENSE.....	17

**PathoSEEK™ Analysis Quick Reference Tables:**
FLOWER

Table 1: Flower Samples ONLY – No Decontamination Step

PathoSEEK™ Assay	Cq Value (High CFU count)	Fluor	Negative Control (Cq)	CFU threshold (CFU/g) (This may vary by state)
<i>Aspergillus</i>	≤ 40	FAM	> 40	Presence/Absence
<i>E. coli</i>	≤ 40	FAM	> 40	Presence/Absence
STEC <i>E. coli</i>	≤ 40	FAM	> 40	Presence/Absence
Salmonella	≤ 40	ROX*	> 40	Presence/Absence
<i>Pseudomonas aeruginosa</i>	≤ 40	FAM	> 40	Presence/Absence
<i>Staphylococcus aureus</i>	≤ 40	FAM	> 40	Presence/Absence
Total Aerobic Count	≤ 20.7	FAM	> 30	100,000 (10 <sup>5</sup> )
Total Coliform (Entero/Col multiplex)	≤ 30.1	FAM	> 40	1,000 (10 <sup>3</sup> )
Total Enterobacteriaceae (Entero/Col multiplex)	≤ 28.1	ROX*	> 40	1,000 (10 <sup>3</sup> )
Total Yeast and Mold	≤ 24.2	FAM	> 40	10,000 (10 <sup>4</sup> )
BTGN (Singleplex)	≤ 28.9	FAM	>40	1,000 (10 <sup>3</sup> )
Total Coliform (BTGN/Coliform Multiplex)	≤ 30.9	ROX	>40	1,000 (10 <sup>3</sup> )
BTGN (BTGN/Coliform Multiplex)	≤ 25.5	FAM	>40	1,000 (10 <sup>3</sup> )
Internal Control*	≤35	HEX	*Internal control verifies the presence or absence of plant DNA	
Assay Positive Control	≤35	FAM/ROX		
Grim Reefer Positive Control	≤35	Cy5		

\*If running singleplex Salmonella or Entero assays, they will be detected on the FAM channel

NON – FLOWER Matrices

Table 2: All Concentrates, MIP Samples – No Decontamination Step (Except gummy, see table 3)

**NOTE:** Grim Reefer does not work in the presence of ALL non flower matrices. See SenSATIVax for MIP SOPs for matrices that have been successfully run with grim reefer

PathoSEEK™ Assay	Cq Value (High CFU count)	Fluor	Negative Control (Cq)	CFU threshold (CFU/g) (This may vary by state)
<i>Aspergillus</i>	≤ 40	FAM	> 40	Presence/Absence
<i>E. coli</i>	≤ 40	FAM	> 40	Presence/Absence
STEC <i>E. coli</i>	≤ 40	FAM	> 40	Presence/Absence
Salmonella	≤ 40	ROX*	> 40	Presence/Absence
<i>Pseudomonas aeruginosa</i>	≤ 40	FAM	> 40	Presence/Absence
<i>Staphylococcus aureus</i>	≤ 40	FAM	> 40	Presence/Absence
Total Aerobic Count	≤ 25.2	FAM	> 35	10,000 (10 <sup>4</sup> )
Total Coliform (Entero/Col multiplex)	≤ 34.7	FAM	> 40	100 (10 <sup>2</sup> )
Total Enterobacteriaceae (Entero/Col multiplex)	≤ 31.8	ROX*	> 40	100 (10 <sup>2</sup> )

PathoSEEK Analysis\_Document\_Agilent\_NODecon\_GR\_v9.1\_DEC2019

Total Yeast and Mold	≤ 37.4	FAM	> 40	1000 (10 <sup>3</sup> )
BTGN (Singleplex)	≤ 33.9	FAM	>40	100 (10 <sup>2</sup> )
Total Coliform (BTGN/Coliform Multiplex)	≤ 35.5	ROX	>40	100 (10 <sup>2</sup> )
BTGN (BTGN/Coliform Multiplex)	≤ 30.9	FAM	>40	100 (10 <sup>2</sup> )
Internal Control*	≤40	HEX	*Internal control verifies the presence or absence of spiked plant positive control (SCCG)	
Assay Positive Controls	≤35	FAM/ROX		
Grim Reefer Positive Control	≤35	Cy5		

\* If running singleplex Salmonella or Entero assays, they will be detected on the FAM channel

### GUMMY

Table 3: Gummy – No Decontamination Step

PathoSEEK™ Assay	Cq Value (High CFU count)	Fluor	Negative Control (Cq)	CFU threshold (CFU/g) (This may vary by state)
<i>Aspergillus</i>	≤ 40	FAM	> 40	Presence/Absence
<i>E. coli</i>	≤ 40	FAM	> 40	Presence/Absence
STEC <i>E. coli</i>	≤ 40	FAM	> 40	Presence/Absence
Salmonella	≤ 40	ROX*	> 40	Presence/Absence
<i>Pseudomonas aeruginosa</i>	≤ 40	FAM	> 40	Presence/Absence
<i>Staphylococcus aureus</i>	≤ 40	FAM	> 40	Presence/Absence
Total Aerobic Count	≤ 27.8	FAM	> 35	10,000 (10 <sup>4</sup> )
Total Coliform (Entero/Col multiplex)	≤ 40	FAM	>40	100 (10 <sup>2</sup> )
Total Enterobacteriaceae (Entero/Col multiplex)	≤ 34.9	ROX*	> 40	100 (10 <sup>2</sup> )
Total Yeast and Mold	≤ 38.1	FAM	> 40	1000 (10 <sup>3</sup> )
BTGN (Singleplex)	≤ 33.7	FAM	>40	100 (10 <sup>2</sup> )
Total Coliform (BTGN/Coliform Multiplex)	≤ 31.7	ROX	>40	100 (10 <sup>2</sup> )
BTGN (BTGN/Coliform Multiplex)	≤ 28.3	FAM	>40	100 (10 <sup>2</sup> )
Internal Control*	≤40	HEX	*Internal control verifies the presence or absence of spiked plant positive control (SCCG)	
Assay Positive Controls	≤35	FAM/ROX		


\* If running singleplex Salmonella or Entero assays, they will be detected on the FAM channel

## Detailed Assay Data Analysis


### 1. Presence / Absence Multiplex Assay: *E. coli* & Salmonella

**NOTE:** If running Salmonella as a singleplex assay, Salmonella will be detected on the FAM channel

- 1.1. Open the Data Analysis window when the run is complete.
- 1.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display
  - Amplification plots will be available for viewing
  - The Cq values will appear to the right in the table
- 1.3. To analyze the results
  - Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

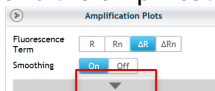

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

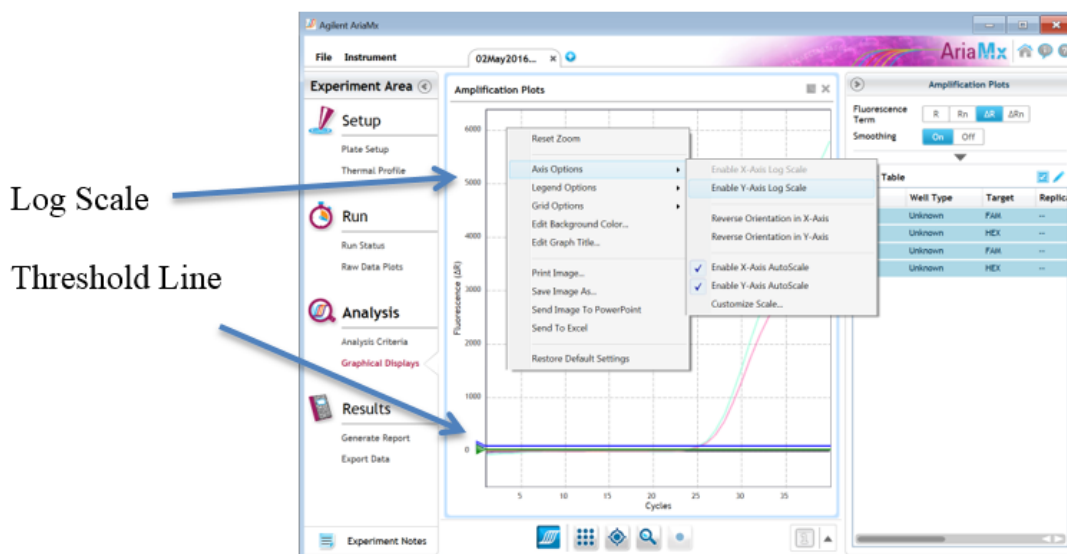

- Controls
  - Assay specific Positive Control, on the FAM and ROX Fluorophores, should have a Cq value  $\leq 35$ .
    - Visually confirm with the curve on the graph.
  - Assay specific Negative Controls, on the FAM and ROX Fluorophores, should have no Cq value.
    - Visually confirm with the curve on the graph.
- Unknown *E. coli* Target (**FAM Fluorophore detects *E. coli***)
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph
  - A “presence” result for the unknown *E. coli* target.
    - Any Cq value for the FAM Fluorophore  $\leq 40$ .
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a presence result occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
  - An “absence” result for the unknown *E. coli* target.
    - No Cq value for the FAM Fluorophore.
    - Visually confirm no curve on the graph.

- **Unknown Salmonella Target (ROX Fluorophore detects Salmonella in multiplex assay, FAM detects Salmonella in singleplex assay)**
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph
  - A “presence” or failing result for the unknown Salmonella target.
    - Any Cq value for the ROX Fluorophore  $\leq 40$ .
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a presence result occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
  - An “absence” or passing result for the unknown Salmonella target.
    - No Cq value for the ROX Fluorophore.
    - Visually confirm no curve on the graph.

## 2. Presence / Absence Singleplex and Multiplex Assay: Aspergillus (includes species specific and multiplex assays)


2.1. Open the Data Analysis window when the run is complete.

2.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

2.3. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

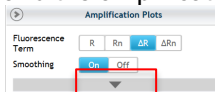

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

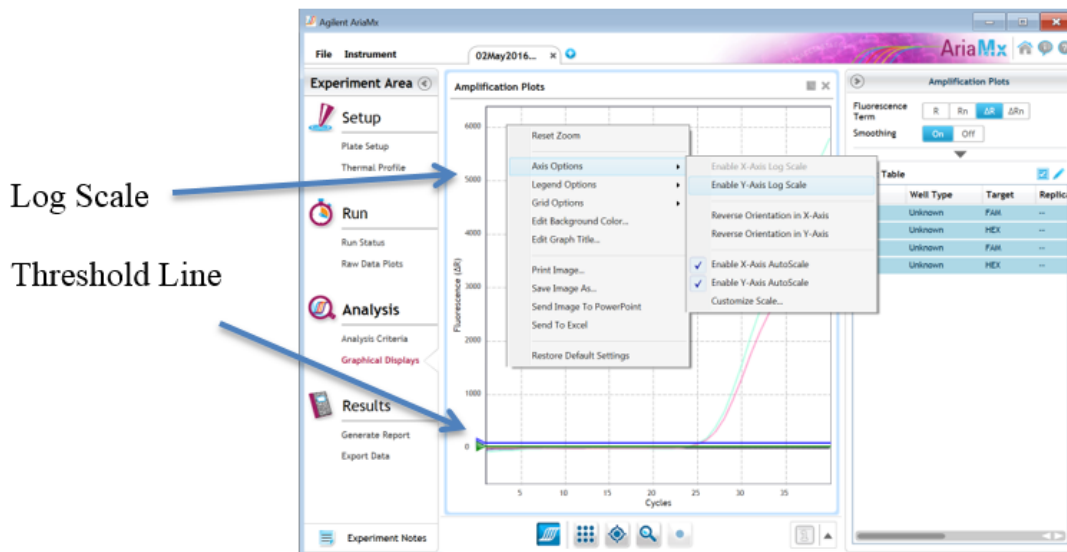

### Controls

- Assay specific Positive Control, on the FAM Fluorophore, has a Cq value  $\leq 35$ .
  - Visually confirm with the curve on the graph.
- Assay specific Negative Control, on the FAM Fluorophore, has no Cq value.
  - Visually confirm with the curve on the graph.
- Grim Reefer Positive Control on the Cy5 Channel, has a Cq value  $\leq 35$ .
- Unknown Aspergillus Target (**FAM Fluorophore detects all 4 Aspergillus species in multiplex assay**)
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph
  - A “presence” or failing result for the unknown Aspergillus target(s).
    - Any Cq value for the FAM Fluorophore  $\leq 40$ .
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a presence result occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
  - An “absence” or passing result for the unknown Aspergillus target.
    - No Cq value for the FAM Fluorophore.
    - Visually confirm no curve on the graph.

### 3. Presence / Absence Multiplex Assay: *Pseudomonas aeruginosa*/Staphylococcus aureus


Open the Data Analysis window when the run is complete.

#### 3.1. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

#### 3.2. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

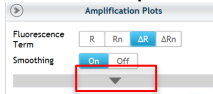

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

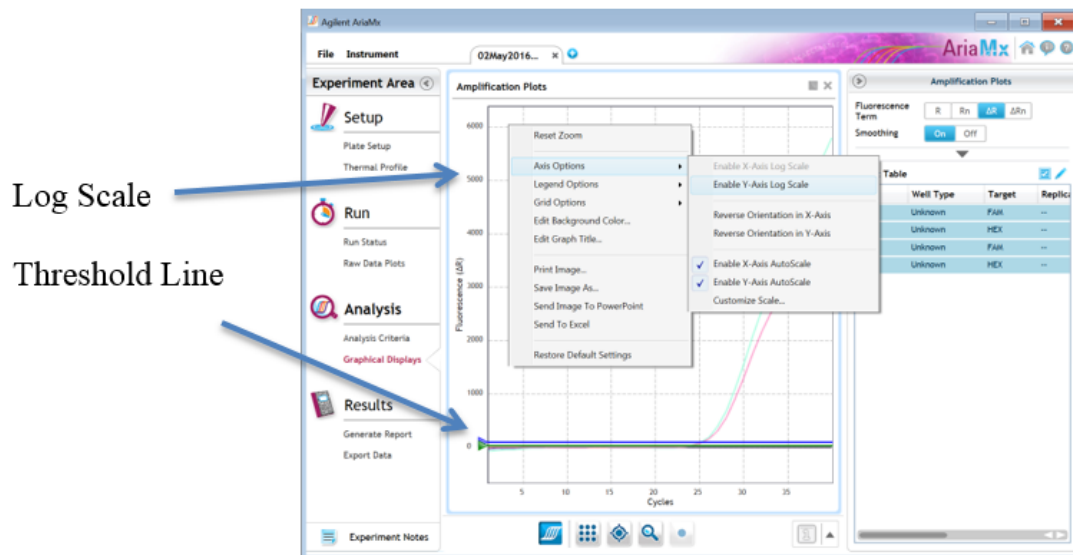

- Controls
  - Assay specific Positive Control, on the FAM Fluorophore, should have a Cq value  $\leq 40$ .
    - Visually confirm with the curve on the graph.
  - Assay specific Negative Control, on the FAM Fluorophore, should have no Cq value.
    - Visually confirm with the curve on the graph.
- Unknown Pseudomonas or Staphylococcus Target (**FAM Fluorophore detects both pathogens**)
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph
  - A “presence” result for the unknown Pseudomonas or Staphylococcus target.
    - Any Cq value for the FAM Fluorophore  $\leq 40$ .
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a presence result occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
  - An “absence” result for the unknown Pseudomonas or Staphylococcus target.
    - No Cq value for the FAM Fluorophore.
    - Visually confirm no curve on the graph.)



#### 4. Threshold Assay: Total Aerobic Count


4.1. Open the Data Analysis window when the run is complete.

4.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

4.3. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

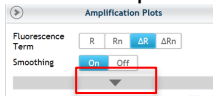

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

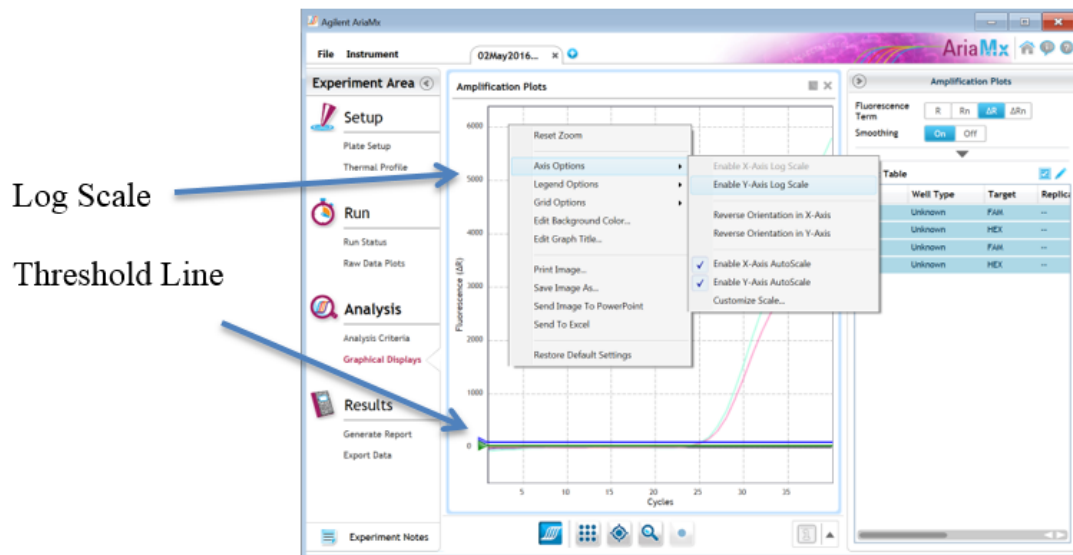

- Controls
  - Assay specific Positive Control, on the FAM Fluorophore, has a Cq value  $\leq 35$ 
    - Visually confirm with the curve on the graph.
  - Assay specific Negative Control, on the FAM Fluorophore, has a Cq value of  $> 30$  or no Cq value.
    - Visually confirm with the curve on the graph.
- Unknown Aerobic Count Target (**FAM Fluorophore detects Total Aerobic Count Bacteria**)
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph.
  - A high CFU count result for the unknown TAC target.
    - **Passing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
  - A low CFU count result for the unknown TAC target.
    - **Failing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph.



## 5. Threshold Assay: Yeast & Mold


5.1. Open the Data Analysis window when the run is complete.

5.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

5.3. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

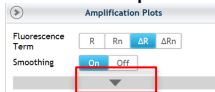

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

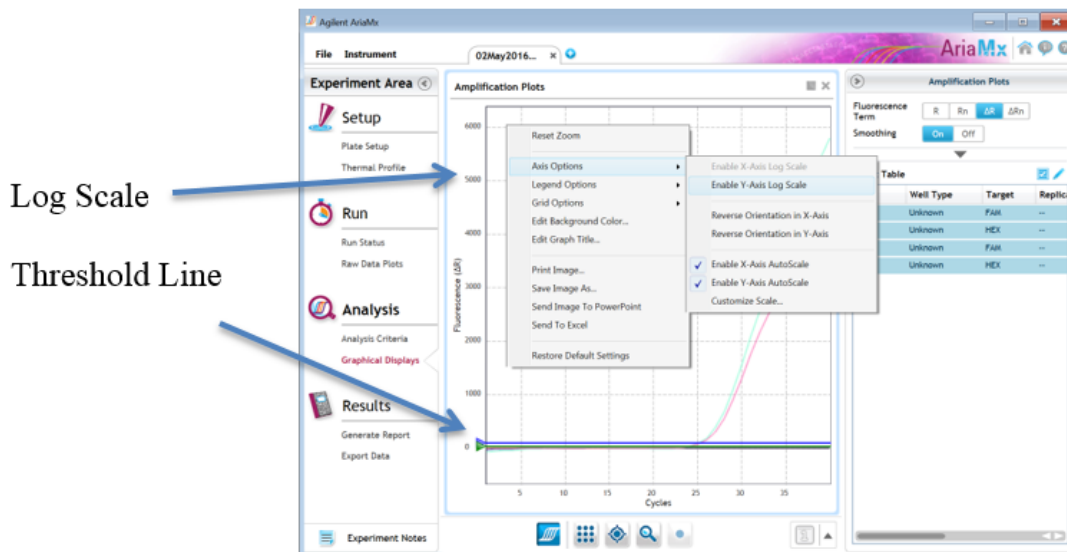

- Controls
  - Assay specific Positive Control, on the FAM Fluorophore, has a Cq value  $\leq 35$ 
    - Visually confirm with the curve on the graph.
  - Assay specific Negative Control, on the FAM Fluorophore, has a Cq value of  $> 35$  or no Cq value.
    - Visually confirm with the curve on the graph.
- Unknown Yeast and Mold Target (**FAM Fluorophore detects Total Yeast & Mold**)
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph.
  - A high CFU count result for the unknown Y&M target.
    - **Passing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
  - A low CFU count result for the unknown Y&M target.
    - **Failing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph.

## 6. Threshold Assay: BTGN – Bile Tolerant Gram-Negative Bacteria


6.1. Open the Data Analysis window when the run is complete.

6.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

6.3. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

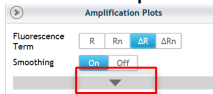

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

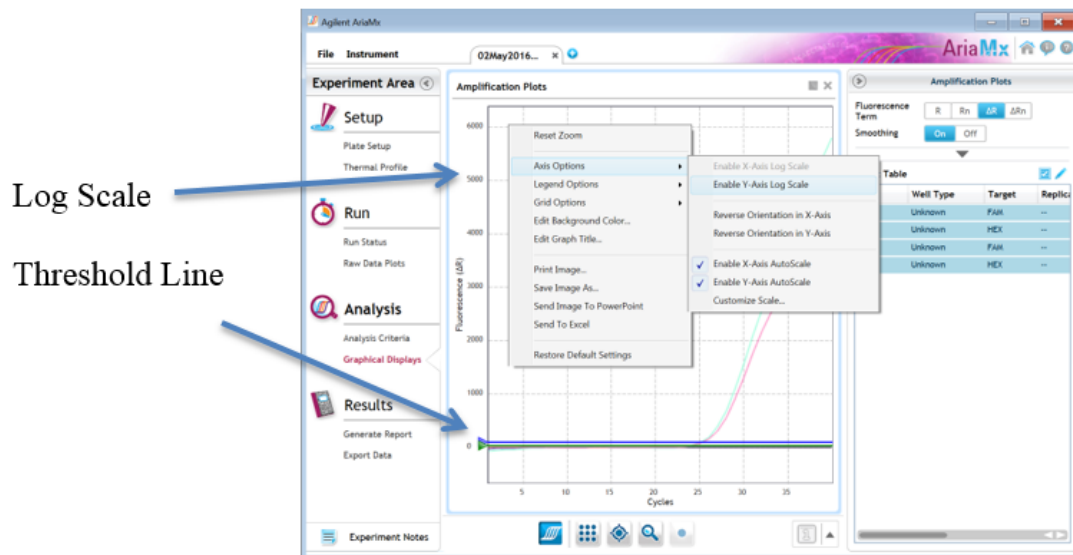

### • Controls

- Assay specific Positive Control, on the FAM Fluorophore, has a Cq value  $\leq 35$ 
  - Visually confirm with the curve on the graph.
- Assay specific Negative Control, on the FAM Fluorophore, has a Cq value of  $> 35$  or no Cq value.
  - Visually confirm with the curve on the graph.

### • Unknown BTGN Target (**FAM Fluorophore detects BTGN**)

- Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
  - Visually confirm with the curve on the graph.
- A high CFU count result for the unknown BTGN target.
  - **Passing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
  - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more details.)
- A low CFU count result for the unknown BTGN target.
  - **Failing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
  - Visually confirm with the curve on the graph.

## 7. Threshold Multiplex Assay: Total Coliform and Entero

**NOTE:** If running Entero as a singleplex assay it will be detected on the FAM channel

7.1. Open the Data Analysis window when the run is complete.

7.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

7.3. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle



- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

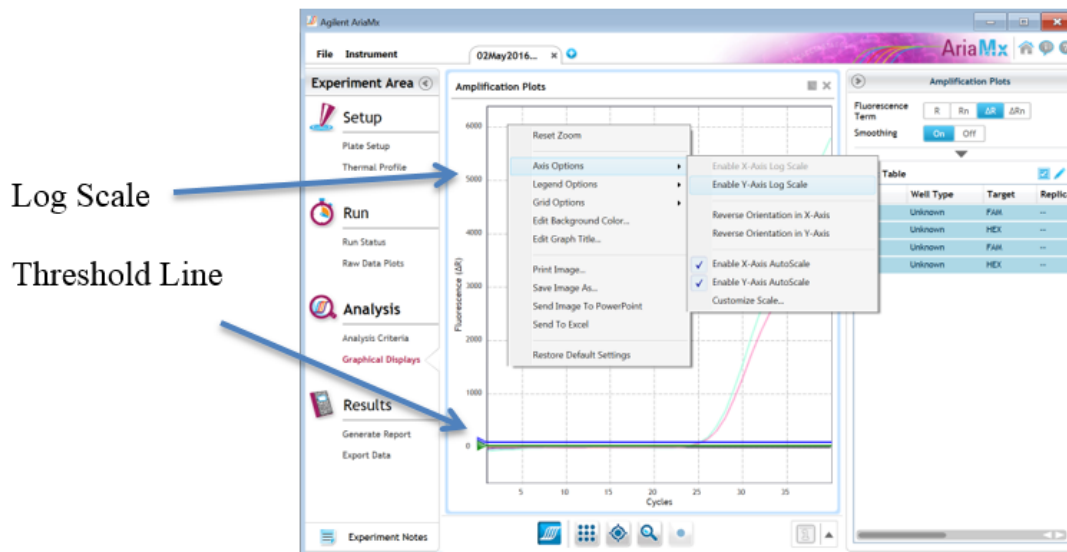

### • Controls

- Assay specific Positive Control, on the FAM and ROX fluorophores, have Cq values  $\leq 35$ 
  - Visually confirm with the curve on the graph.
- Assay specific Negative Control, on the FAM and ROX fluorophores, have Cq values  $> 35$  or no Cq value.
  - Visually confirm with the curve on the graph.

### • Unknown Coliform Target (**FAM Fluorophore detects Total Coliform**)

- Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
  - Visually confirm with the curve on the graph.
- A high CFU count result for the unknown coliform target.
  - **Passing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
  - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more detail.)
- A low CFU count result for the unknown coliform target.
  - **Failing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
  - Visually confirm with the curve on the graph.

- Unknown Entero Target (ROX Fluorophore detects Entero in the multiplex assay, FAM Fluorophore detects Entero in the singleplex assay)
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for non-flower matrices.
    - Visually confirm with the curve on the graph.
  - A high CFU count result for the unknown Entero target.
    - Passing Sample Result:** Check Cq Value on the ROX Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. -See troubleshooting guide below for more detail.)
  - A low CFU count result for the unknown Entero target.
    - Failing Sample Result:** Check Cq Value on the ROX Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph.

## 8. Threshold Multiplex Assay: BTGN and Total Coliform

**NOTE:** If running Coliform or BTGN a singleplex assay it will be detected on the FAM channel


8.1. Open the Data Analysis window when the run is complete.

8.2. Highlight the wells of interest in the Analysis Criteria under Analysis, then select Graphical Display


- Amplification plots will be available for viewing
- The Cq values will appear to the right in the table

8.3. To analyze the results

- Start by turning the graph to Log Scale with a right click on the chart, select Axis options, enable y-axis log scale. Expand the amplification plots settings by clicking on the triangle

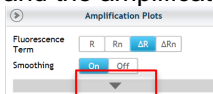

- Manually adjust thresholds to 100 RFU for the HEX, FAM, ROX and Cy5 fluorophores.

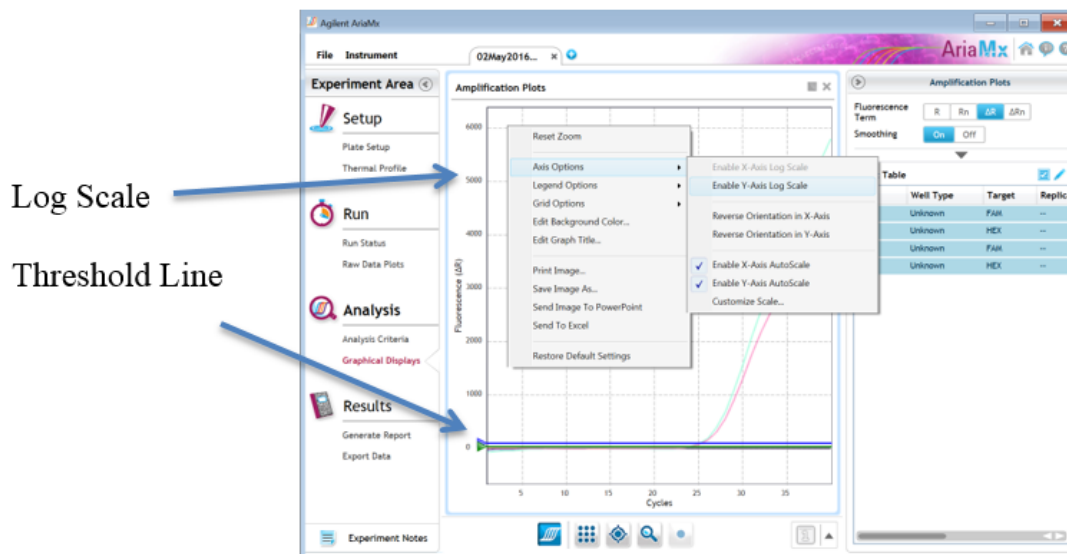

- Controls
  - Assay specific Positive Control, on the FAM and ROX fluorophores, have Cq values  $\leq 35$ 
    - Visually confirm with the curve on the graph.
  - Assay specific Negative Control, on the FAM and ROX fluorophores, have Cq values  $> 35$  or no Cq value.

- Visually confirm with the curve on the graph.
- **Unknown Coliform Target (ROX Fluorophore detects Total Coliform)**
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for all other matrices.
    - Visually confirm with the curve on the graph.
  - A high CFU count result for the unknown coliform target.
    - **Passing Sample Result:** Check Cq Value on the ROX Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. (See troubleshooting guide below for more detail.)
  - A low CFU count result for the unknown coliform target.
    - **Failing Sample Result:** Check Cq Value on the ROX Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph.
- **Unknown BTGN Target (FAM Fluorophore detects BTGN in the multiplex assay)**
  - Internal Control, on the HEX Fluorophore, has a Cq value  $\leq 35$  for flower samples,  $\leq 40$  for non-flower matrices.
    - Visually confirm with the curve on the graph.
  - A high CFU count result for the unknown BTGN target.
    - **Passing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph. It is very important to confirm with the amplification curve when a high CFU count occurs. Sometimes the background amplification will give a false positive reading, especially when Cq reading is less than 15. -See troubleshooting guide below for more detail.)
  - A low CFU count result for the unknown BTGN target.
    - **Failing Sample Result:** Check Cq Value on the FAM Fluorophore. See Tables 1-3 for Cq cutoff value depending on matrix being tested.
    - Visually confirm with the curve on the graph.

**Table 4. Cq to CFU Conversion Equation Table**

Matrix	Microbial Test	Cq to CFU/g Conversion Equation
Flower	Total Yeast and Mold	$CFU/g = 10^{[(36.671 - Cq \text{ Value})/3.1194]}$
Flower	Total Aerobic Count	$CFU/g = 10^{[(35.111 - Cq \text{ Value})/2.8883]}$
Flower	Total Coliform (Singleplex and Ent/Col MPX)	$CFU/g = 10^{[(40.073 - Cq \text{ Value})/3.3417]}$
Flower	Total Enterobacteriaceae (Singleplex and Ent/Col MPX)	$CFU/g = 10^{[(41.218 - Cq \text{ Value})/4.3708]}$
Flower	BTGN (singleplex)	$CFU/g = 10^{[(38.665 - Cq \text{ Value})/3.2417]}$
Flower	BTGN (BTGN/Coliform MPX)	$CFU/g = 10^{[(34.908 - Cq \text{ Value})/3.1506]}$
Flower	Coliform (BTGN/Coliform MPX)	$CFU/g = 10^{[(44.665 - Cq \text{ Value})/4.5877]}$
MIP/Extract	Total Yeast and Mold	$CFU/g = 10^{[(54.972 - Cq \text{ Value})/5.8485]}$
MIP/Extract	Total Aerobic Count	$CFU/g = 10^{[(38.076 - Cq \text{ Value})/3.2249]}$
MIP/Extract	Total Coliform (Singleplex and Ent/Col MPX)	$CFU/g = 10^{[(41.935 - Cq \text{ Value})/3.6274]}$
MIP/Extract	Total Enterobacteriaceae (Singleplex and Ent/Col MPX)	$CFU/g = 10^{[(38.407 - Cq \text{ Value})/3.3041]}$
MIP/Extract	BTGN (singleplex)	$CFU/g = 10^{[(40.58 - Cq \text{ Value})/3.3403]}$
MIP/Extract	BTGN (BTGN/Coliform MPX)	$CFU/g = 10^{[(38.623 - Cq \text{ Value})/3.8199]}$
MIP/Extract	Coliform (BTGN/Coliform MPX)	$CFU/g = 10^{[(45.591 - Cq \text{ Value})/5.015]}$
Gummy	Total Yeast and Mold	$CFU/g = 10^{[(52.989 - Cq \text{ Value})/4.9718]}$
Gummy	Total Aerobic Count	$CFU/g = 10^{[(37.235 - Cq \text{ Value})/2.356]}$
Gummy	Total Coliform (Singleplex and Ent/Col MPX)	$CFU/g = 10^{[(52.888 - Cq \text{ Value})/5.9643]}$
Gummy	Total Enterobacteriaceae (Singleplex and Ent/Col MPX)	$CFU/g = 10^{[(44.81 - Cq \text{ Value})/4.9665]}$
Gummy	BTGN (singleplex)	$CFU/g = 10^{[(40.458 - Cq \text{ Value})/3.4035]}$
Gummy	BTGN (BTGN/Coliform MPX)	$CFU/g = 10^{[(34.483 - Cq \text{ Value})/3.0631]}$
Gummy	Coliform (BTGN/Coliform MPX)	$CFU/g = 10^{[(39.366 - Cq \text{ Value})/3.7968]}$


Please Contact [support@medicinalgenomics.com](mailto:support@medicinalgenomics.com) for an easy to use conversion spreadsheet



## 9. Export the Data


### 9.1. Exporting the Cq values into an Excel spreadsheet.

- To export the Cq values to an Excel spreadsheet, right-click on the chart on the bottom right of the screen.
- Choose Send To Excel...

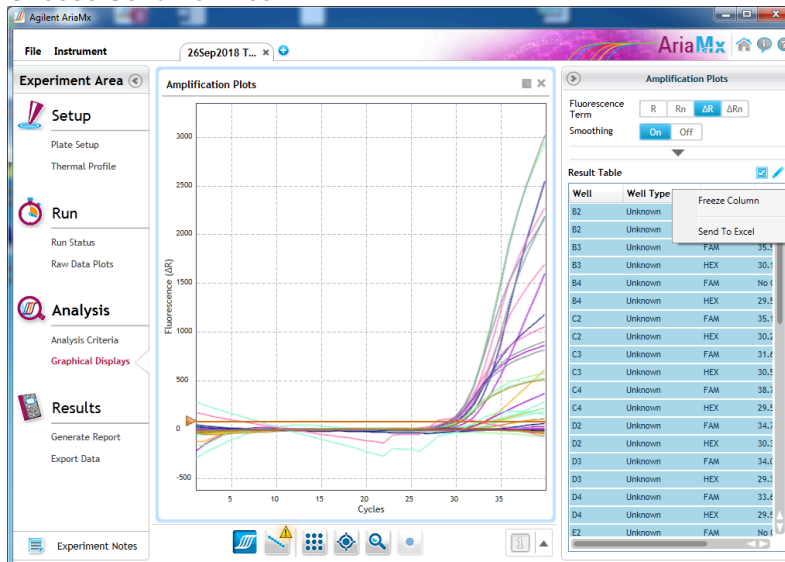


### 9.2. Saving a visual of the graph

- To save a picture of the graph, navigate to the Results section of the software and select Generate Report
- Select PDF or Powerpoint and select Generate Report to export pictures

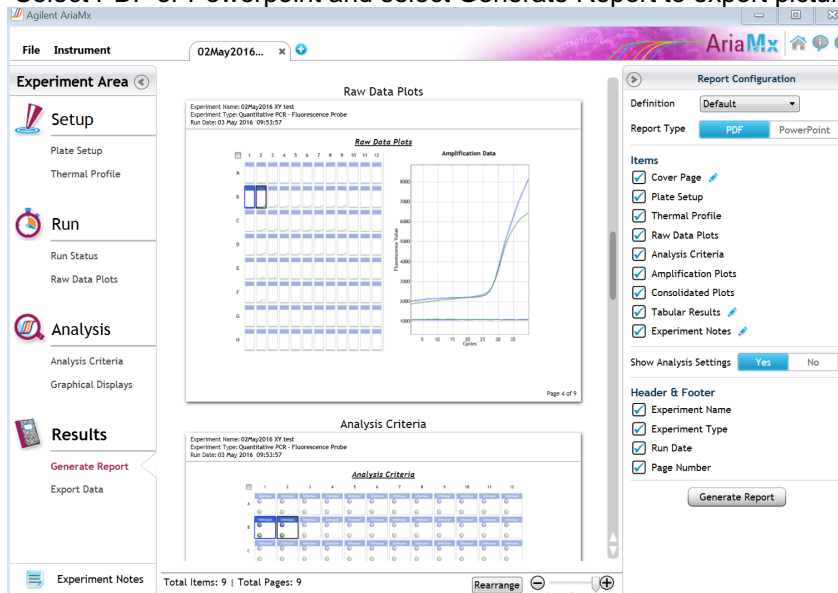



## Troubleshooting Guide

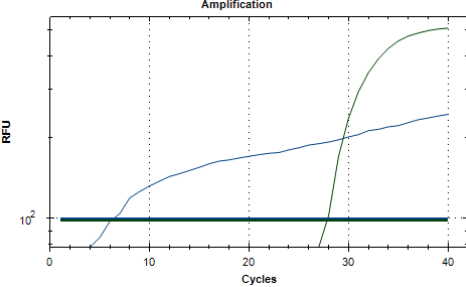Symptom	Reason	Solution
Internal control (SCCG Primer) failure	Extraction Failure	Repeat SenSATIVAx™ and PathoSEEK™ by following the protocol.
	Residual ethanol in elution	Ethanol is an inhibitor to PCR. Return to the SenSATIVAx™ protocol and repeat all steps.
	Mix up in Reaction Setup	Repeat the qPCR by following the protocol.
	Missing Fluorophore on plate set up	In the Data Analysis window click on View/Edit Plate Setup from the Settings drop down. All wells should have both FAM and HEX. Once completed and window is closed the analysis should automatically update.
Internal Control (SCCG) Positive result on positive or negative control samples or samples that do not contain plant DNA <b>No Cq or a Cq of more than 35 is acceptable. Any Cq lower than 35 constitutes a rerun</b>	Plant DNA contamination in a reagent	Troubleshoot which reagent was contaminated; use new reagents, thoroughly clean all pipettes and bench areas with 10% bleach solution.
	qPCR bench too close to extraction area	Designate separate benches, pipettes etc. for extractions and qPCR setup
Positive Negative Control	Small Cq value <15	Visually confirm that there is an amplification curve. If not, this is low level background and is to be expected.
	Carry over	Repeat the qPCR by following the protocol.
	Insufficient pre-setup bleaching	Wipe down the lab workspace and all equipment with 10% Bleach. Repeat qPCR.
Negative Positive Control	Mix up in Reaction Setup	Repeat the qPCR by following the protocol.
Total run failure	Excessive vortex of the qPCR Master Mix	Repeat the qPCR by following the protocol.
Background Amplification 	Unclear	This is usually seen with a very low Cq reading (<15), the curve is usually missing the exponential growth phase, but rather a gradual increase of fluorescent signal. This is usually a negative result, but should be repeated.
No Cy5 Signal	Did not use Grim Reefer Positive control	Add Grim Reefer Positive Control to sample during SenSATIVAx DNA Extraction

## **Glossary and Definitions**

**Deoxyribonucleic acid (DNA)** is a molecule that encodes the genetic instructions used in the development and functioning of all known living organisms.

**Polymerase Chain Reaction (PCR)** is a technology in molecular biology used to amplify a single copy or a few copies of a piece of DNA across several orders of magnitude, generating thousands to millions of copies of a particular DNA sequence.

A **fluorophore** is a fluorescent chemical compound that can re-emit light upon light excitation.

The **Negative Controls** are the reactions where no Cq is expected. It helps to ensure that all Assay specific reactions are clean of contaminants.

The assay specific **Positive Controls** are the reactions where a Cq is expected. It helps ensure that all Assay specific reactions are working correctly. The Assay specific Positive Control is targeting the pathogen using the FAM Fluorophore.

The **Internal Control** is added to every sample reaction where a Cq is expected. It ensures the effectiveness and efficiency of each reaction. The internal control targets plant DNA, or more specifically, a Single Copy Control Gene (SCCG), using the HEX Fluorophore.

**MIP** is short for Marijuana Infused Product. A MIP is cannabis plant material or concentrate mixed into a consumable.

### **DISCLAIMER**

This test was developed, and its performance characteristics determined by Medicinal Genomics Company, for laboratory use. Any deviations from this protocol are not supported by MGC.

The results may vary based on laboratory conditions. Altitude and humidity are among factors known to affect the growth of bacterial and fungal species. All thresholds were determined based on the results using the BIO-RAD CFX96 Touch™ Real-Time PCR Detection System. It is recommended that thresholds be calibrated for each specific laboratory setting.

### **LIMITED USE LABEL LICENSE**

This product is covered by at least one or more claims of US patents applications, which are exclusively licensed to Medicinal Genomics Corporation. This product is sold strictly for the use of the buyer, and the buyer is not authorized to transfer this product [or any materials made using this product] to any third party.

© 2019 Medicinal Genomics Corporation. All rights reserved.

\* All Trademarks are property of their respective owners.